



BDL-SMN

BDL-SMN Picosecond / CW Diode Laser Family

Free-beam output or single-mode fibre coupling

Beam-profile correction optics

Wavelengths 375 nm, 405 nm, 445 nm, 473 nm, 488 nm, 515 nm, 640 nm, 685 nm, 785 nm

Pulsed and CW operation

Pulse width down to 40 ps

Repetition rate 20-50-80 MHz

Low skew trigger output

Cooled laser diode

Internal power regulation loop

Linear response to power control signal

Fast on / off / multiplexing capability

Synchronisation input

Complete electronics integrated in laser housing

Simple +12V wall-mounted power supply


Luminescence lifetime experiments

Laser scanning microscopy

Fluorescence correlation

Time-correlated single photon

counting experiments

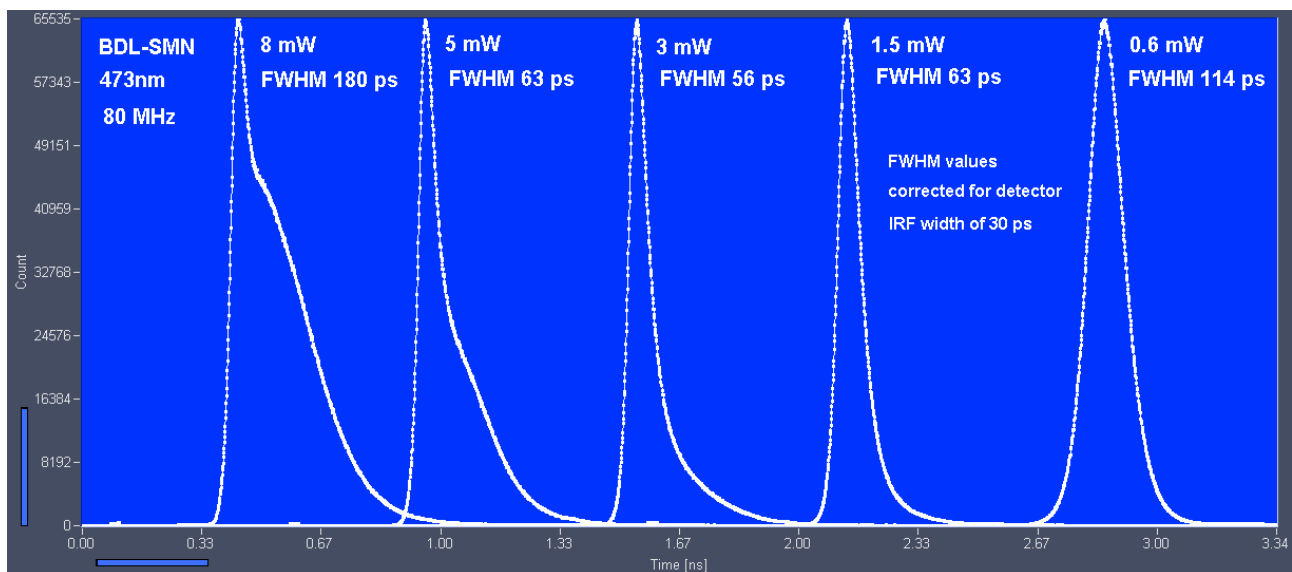

Designed and manufactured by



Becker & Hickl GmbH
Nahmitzer Damm 30
12277 Berlin, Berlin
Tel. +49 / 30 / 787 56 32
Fax. +49 / 30 / 787 57 34
email: info@becker-hickl.com
www.becker-hickl.com



LASOS Lasertechnik GmbH
Carl-Zeiss-Promenade 10
07745 Jena, Germany
Tel. +49 3641 2944-0
Fax +49 3641 2944-17
info@lasos.com
www.lasos.com



BDL-SMN

Optical

Repetition Rate
Wavelength, nm
Pulse width (FWHM, at medium power)
Pulse width (FWHM, at maximum power)
Peak Power
Power control range
(Average CW equivalent power,
adjustable via external power control signal)


Diameter of laser beam
Polarisation
Fibre coupling
Coupling efficiency into single-mode fibre, typically
Stability of Repetition Rate
Pulse-to Pulse Jitter
Reaction time to 'Laser on' signal (pulsed mode)
Reaction time to 'Laser on' signal (CW mode)
Power and pulse shape stabilisation after switch-on

20-50-80 MHz, or CW operation
375, 405, 445, 473, 488, 515, 640, 685, 785, other on request
40 to 90 ps ²⁾
200 to 300 ps ²⁾
40 to 500 mW ¹⁾
20 MHz: 0 to 0.6 mW 0 to 2 mW ²⁾
50 MHz: 0 to 1.5 mW 0 to 5 mW ²⁾
80 MHz: 0 to 2.4 mW 0 to 8 mW ²⁾
CW mode: 0 to 20 mW 0 to 50 mW ²⁾
0.7 mm, TEM₀₀ mode
horizontal
Kineflex system of Qioptiq
60%
± 100 ppm
< 20 ps
3 µs
3 µs
2 min ⁵⁾

Trigger Output

Pulse Amplitude
Pulse Width
Output Impedance
Connector
Delay from Trigger to Optical Pulse
Jitter between Trigger and Optical Pulse

1 V (peak) into 50 Ω
1 ns
50 Ω
SMA
< 1 ns
< 10 ps

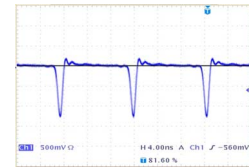

Synchronisation Input

Amplitude
Duty cycle
Frequency
Switching from internal clock to snc input

+3.3 to +5V into 50 Ω
10 to 30 %. DC equivalent must be < 2.5V
20 to 80 MHz
Automatic, by average voltage at sync input connector

Control Inputs

Frequency 20 MHz
Frequency 50 MHz
Frequency 80 MHz
CW operation
Laser ON / Off
External Power Control

TTL / CMOS high ³⁾
TTL / CMOS high ³⁾
TTL / CMOS high ³⁾
TTL / CMOS high ³⁾
TTL / CMOS low ³⁾
analog input, 0 to + 10V

Power Supply

Power Supply Voltage
Power Supply Current
Power Adapter

+ 9 V to +12 V
300 mA to 1.5 A ⁴⁾
AC-DC power adapter, with key switch and control box in cable

Mechanical Data

Dimensions
Mounting Thread

160 mm x 90 mm x 60 mm
two M6 holes

Maximum Values

Power Supply Voltage
Voltage at Digital Control Inputs
Voltage at Ext. Bias Input
Ambient Temperature

0 V to +15 V
-2 V to +7 V
-12 V to + 12 V
0 °C to 40 °C ⁵⁾

- 1) Typical values, sample tested. Depends on pulse width and selected power.
- 2) Depends on wavelength version.
- 3) All inputs have 10 kΩ pull-up resistors. Open input is equivalent to logic 'high'.
- 4) Dependent on ambient temperature. Cooling current changes due to temperature regulation of laser diode
- 5) Operation below 13 °C may result in extended warm-up time.



Caution: Class 3B laser product. Avoid direct eye exposure. Light emitted by the device may be harmful to the human eye. Please obey laser safety rules when operating the devices. Complies with US federal laser product performance standards.

International Sales Representatives




US:
Boston Electronics Corp
tcspc@boselec.com
www.boselec.com



UK:
Photonic Solutions PLC
sales@psplc.com
www.psplc.com



Japan:
Tokyo Instruments Inc.
sales@tokyoinst.co.jp
www.tokyoinst.co.jp

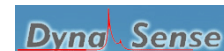

China:
DynaSense Photonics Co. Ltd.
info@dyna-sense.com
www.dyna-sense.com